

ThP 099 NanoMEA Chip Platform for Proteomics

When: Thursday, June 9th; 10:30am – 2:30pm CT

Poster presenter: Robert Maxwell, Ph.D., Senior Scientist at Newomics Inc.

Introduction

Proteomics of small volumes of biological samples down to single cells has progressed rapidly. However, the sensitivity and reproducibility of ultralow flow LC-MS for proteomics has remained a challenge. We developed a silicon-based, plug and play, nanoflow Multinozzle Emitter Array (nanoMEA) chip to directly address this challenge. NanoMEA chip monolithically integrates our M3 multi-nozzle emitter and an on-chip LC column, thereby reducing the dead volume and simplifying the plumbing and connection for ultraflow LC-MS. Herein, we benchmark the performance of nanoMEA chip for small-volume top-down and bottom-up proteomics.

Methods

NanoMEA chips were designed using the L-Edit software (v15, Tanner Research Inc.) and fabricated as we published before [Anal.Chem.2011,83,6082–6089]. The fabricated devices were examined by optical microscopy using a Reichert-Jung Polylite 88 microscope (Reichert Microscope Services) and by scanning electron microscopy using a JEOL 6340F FEG-SEM (JEOL Ltd.). On-chip columns were packed with C18 and C4 beads for bottom-up and top-down proteomics applications, respectively. Nanoflow LC-MS and proteomic analysis were performed using a Newomics MnESI ion source interfaced to a Thermo Orbitrap Eclipse Tribrid with a Vanquish Neo UPLC, or a TIMS-PASEF TOF equipped with a nanoElute. Standard proteins and their tryptic digests as well as HeLa digests were used for testing the performance of our nanoMEA chips.

Preliminary Data

We developed a prototype MEA chip with 10 μm single and double emitters and an on-chip LC column. The columns have 75 μm and 150 μm id with a length ranging from 5 cm to 20 cm. The chips were tested with a variety of chromatography packing beads, including 1.9 μm C18 packing for RPLC and 3 μm C4 packing for intact protein analysis. We achieved 3 second peak widths using C18 packing for peptide 722 (m/z) from BSA digests. A comparison was made for a nanoMEA chip with a chip based internal column: 75 μm X 15cm and 1.9 μm packing. This prototype chip was developed with one nozzle and with two nozzles. We show that an 50-80% increase in sensitivity when we used two nozzles rather than 1 nozzle. As such, two nozzles splits the 400 nl/min flow rate for each 10 μm nozzle to 200 nl/min, thereby increasing the sensitivity through increase in ESI efficiency. We compared the nanoMEA chip prototype to traditional ESI in details. A fused silica column with the same column dimensions and packing was produced in house and connected to a 10 μm silica Picotip emitter tip. We found that the single nozzle nanoMEA Chip had a 25% lower peak width and 50% higher sensitivity for BSA digests and a double nozzle MEA chip had nearly 2-fold higher overall sensitivity. Proteomic analysis for the 2 nozzle nanoMEA chip yielded 3 fold higher peptide counts for lower amounts of Hela digest when compared to the traditional ESI.

Novel Aspect

First demonstration of a monolithic silicon microfluidic chip that integrates a LC column with multi-nozzle emitter for ultralow flow proteomics.